

INTEGRATING NEPHELOMETER MODEL 3563:RP-3563

STANDARD SERVICE PROCEDURES

Standard maintenance and calibration service for the Integrating Nephelometer Model 3563 consists of these procedures:

- 1) Review written instructions or explanation of problem from customer
- 2) Initial inspection
- 3) Update manual, software and firmware as necessary
- 4) Clean instrument and optics
- 5) Inspect and replace as needed: filters, tubing, and O-rings
- 6) Leak check
- 7) Align optics, apertures and shutters
- 8) Calibrate humidity and temperature sensors
- 9) Calibrate unit with filtered air and CO₂
- 10) Final electrical checkout and performance verification

Standard service includes minor repairs and replacement parts. Major repairs and replacement parts will be charged for parts and labor. TSI will contact the customer for approval before making any major repairs.

Turnaround time for standard service is approximately two weeks after receipt of the instrument. TSI service work is backed by a 90 day service warranty.



UNDERSTANDING, ACCELERATED

TSI Incorporated – Visit our website www.tsi.com for more information.

USA	Tel: +1 800 874 2811	India	Tel: +91 80 67877200
UK	Tel: +44 149 4 459200	China	Tel: +86 10 8251 6588
France	Tel: +33 4 91 11 87 64	Singapore	Tel: +65 6595 6388
Germany	Tel: +49 241 523030		