

ThermalPro will not allow Velocity to be Analyzed

Frequently Asked Question #14

Question:

I have acquired data in ThermalPro. When I try to “Analyze” my data, it will not allow me to analyze “velocity” – It says “velocity over range.” When I analyze the text file and then look at it, only the first ½ of the samples are there, the second half are all zeroes. What is the problem?

Answer:

The DAC driver is corrupted. Re-install the DAC driver using the following method:

1. Go to the Control Panel->System->Hardware->Device Manager.
2. Pick the UEI device from the list.
3. Right mouse click on the UEI device to bring up the “Properties” page.
4. Go to the “Driver” page.
5. Click “Update Driver”
6. Follow the driver installation instruction; select the ThermoPro CD so Windows can search for the driver.



TSI Incorporated - 500 Cardigan Road, Shoreview, MN 55126-3996 USA
USA Tel: +1 800 874 2811 E-mail: fluid@tsi.com Website: www.tsi.com
UK Tel: +44 149 4 459200 E-mail: tsuuk@tsi.com Website: www.tsiinc.co.uk
France Tel: +33 491 95 21 90 E-mail: tsifrance@tsi.com Website: www.tsiinc.fr
Germany Tel: +49 241 523030 E-mail: tsigmbh@tsi.com Website: www.tsiinc.de
Sweden Tel: +46 8 595 13230 E-mail: tsiab@tsi.com Website: www.tsi.se
India Tel: +91 80 41132470 E-mail: tsi-india@tsi.com
China Tel: +86 10 8260 1595 E-mail: tsibeijing@tsi.com

Contact your local TSI Distributor or visit our website www.tsi.com for more detailed specifications.



TRUST. SCIENCE. INNOVATION.