

A component-based system to produce monodisperse, submicrometer particles

Using an Electrostatic Classifier 3082, the model 3740A extracts a narrow monodisperse size range from polydisperse aerosol produced with an atomizer. The resulting monodisperse aerosol will typically have a geometric standard deviation of 1.05 micrometer or better.

Features and Benefits

- Generates a highly monodisperse aerosol in the size range from 0.01 to 1.0 micrometer
- Operator simply selects the operating parameters required to produce a known particle size
- Operation is based on first principles, so no calibration is necessary

Applications

This complete aerosol generation system is useful anywhere a precise, monodisperse, submicrometer aerosol is needed. Specific applications include:

- Calibrating submicrometer counting and sizing instruments (Condensation Particle Counters and Optical Particle Counters) as described in JIS B 9921
- Testing HEPA, ULPA, and other high-efficiency filters as described in IES RP-CC007.1 and RP-CC021.1, and DIN 24183
- Studying nucleation, condensation, and diffusion
- Studying particle charge and electrical mobility
- Enhancing the monodispersity of polystyrene latex (PSL) and other aerosols



Model 3940A Operation Overview

Operation

Compressed air first passes through a Filtered Air Supply, where the air is cleaned and dried. This conditioned air is used to power a Constant Output Atomizer. The atomizer generates an aerosol of polydisperse droplets with a count median diameter of 0.3 micrometer. It works with oils such as DOP and DEHS, most solutions with a volatile solvent, or suspensions such as PSL spheres in water. The aerosol droplets pass through a Diffusion Dryer where any volatile components evaporate off. The remaining particles then pass through Kr-85 bipolar chargers to establish a low-level charge equilibrium.

The Electrostatic Classifier used in the system contains a Differential Mobility Analyzer (DMA). Inside the DMA, particles are subjected to an electric field. The field deflects the particles from their flowpath according to their electrical mobility, which is inversely proportional to particle size. By adjusting voltage levels, particles of a narrow size range are steered into a small exit. The particles leave the classifier as a monodisperse aerosol. Aerosol concentration is changed simply by adjusting valves so that a portion of the aerosol passes through high-efficiency filters. (For more information, see the product information sheet of the Electrostatic Classifier 3082).


Bibliography

Caldow R and J Blesener, A Procedure to Verify the Lower Counting Limit of Optical Particle Counters, *Journal of Parenteral Science and Technology*, July/August 1989. (TSI paper A62)

Kinney PK, DYH Pui, GW Mulholland, and NP Bryner, Use of the Electrostatic Classification Method to Size 0.1µm SRM Particles—A Feasibility Study, *Journal of Research of the National Institute of Standards and Technology* 96(2), March/April 1991. (TSI paper A74)

Liu BYH and DYH Pui, A Submicron Aerosol Standard and the Primary, Absolute Calibration of the Condensation Nuclei Counter, *Journal of Colloid and Interface Science* 47(1), April 1974. (TSI paper A81)

Liu BYH, DYH Pui, and KY Rubow, Performance of HEPA and ULPA Filters, Proceedings of the 1985 annual technical meeting of the Institute of Environmental Sciences. (*TSI paper A47*)

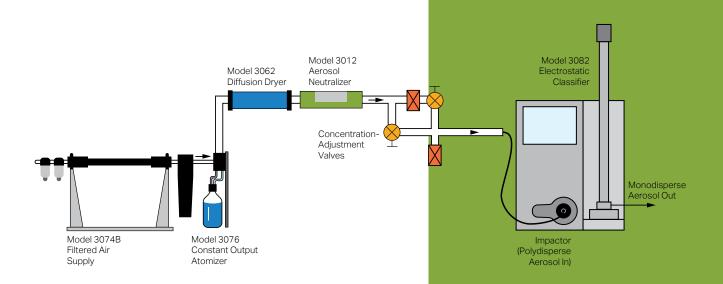

Optional Configuration

The standard model 3940A system uses two neutralizers to ensure that particle charge is fully reduced to a Fuch's equilibrium before aerosol enters the DMA. By itself, the 3077A may not be adequate if the high-concentration aerosol generated by the atomizer is also highly charged.



This 3940A system is shown ready to calibrate a TSI® Electrostatic Classifier 3082. Ask your TSI® representative about additional components needed for your particular application.



Specifications

Submicrometer Monodisperse Aerosol Generation System Model 3940A

Refer to separate product sheets for descriptions and specifications of individual system components.

Equipment Included

3082 Electrostatic Classifier Platform

3081A Long DMA
3077A Aerosol Neutralizer
3012 Aerosol Neutralizer
3074B Filtered Air Supply
3076 Constant Output Atomizer

3062 Diffusion Dryer

Particle Material

Solid or non-volatile liquid

Particle Size Range

0.005 to $1 \, \mu m$

Particle Concentration

Adjustable from 0 to 10⁵ particles/cm³

Charger

Bipolar, Kr-85, 2 millicurie, half-life of 10.4 years*

Aerosol Flow Rate

0.2 to 3.5 L/min

Aerosol Pressure Range at Classifier Outlet

1 ± 0.2 atm

Monodispersity

±5% of mean size for singly charged particles

Compressed-Air Requirements

25 L/min at 35 psi

* TSI is authorized by the United States Nuclear Regulatory Commission to distribute these neutralizers to customers in most countries. If your location is within the United States, no other federal license is required. Check local regulations for your own protection.

Specifications are subject to change without notice.

TSI and the TSI logo are registered trademarks in the United States and may be protected under other country's trademark registrations

To Order

Submicrometer Monodisperse Aerosol Generation System

Specify Description

3940A Submicrometer Monodisperse Aerosol Generation System

Optional Accessories

Specify Description

3001788 Conductive tubing, 1/4-inch inside diameter 3001789 Conductive tubing, 3/8-inch inside diameter

3068B Aerosol Electrometer 3708 Flow Splitter

3750 Condensation Particle Counter*

*Other CPCs may be used



 $\textbf{TSI Incorporated -} \ \textit{Visit our website } \textbf{www.tsi.com} \ \textit{for more information}.$

 USA
 Tel: +1 800 874 2811
 India
 Tel: +91 80 67877200

 UK
 Tel: +44 149 4 459200
 China
 Tel: +86 10 8219 7688

 France
 Tel: +33 1 41 19 21 99
 Singapore
 Tel: +65 6595 6388

 Germany
 Tel: +49 241 523030

P/N 5001619 (A4) Rev F

©2022 TSI Incorporated

Printed in U.S.A.